



Specifications				
Item		CX41		
Optical system		UIS2 (Universal Infinity-corrected) optical system		
Illumination		Built-in transmitted Koehler illuminator 6 V 30 W halogen bulb 100–120 V/220–240 V~ 0.85/0.45 A 50/60 Hz		
Focusing		●Stage height movement by roller guide (rack & pinion) ●Stroke per rotation: 36.8 mm ●Full stroke range: 25 mm ●Upper limit stopper ●Tension adjustment on coarse focus adjustment knob		
Revolving nosepiece		Fixed quintuple nosepiece with inward tilt		
Observation tube	Type	U-CBI30-2, Binocular	U-CTR30-2, Trinocular	U-CTBI, Tilting binocular
	Field number	20	20	18
	Tube inclination	30°	30°	30°—60°
	Interpupillary distance adjustment range	48—75 mm	48—75 mm	48—75 mm
	Light path selector	None	None (Bi 50%, Video/Photo 50% fixed)	None
Stage	Size	188(W) x 134(D) mm		
	Movement range	76 mm X-direction x 50 mm Y-direction		
	Specimen holder	Double slide holder		
	Rubber grip	Equipped as standard		
Condenser (CH3-CD)	Type	Abbe condenser		
	N.A.	1.25 with oil immersion		
	Aperture iris diaphragm	Built-in		
Dimensions & weight		233(W) x 432(H) x 367.5(D) mm, approximately 8.5 kg (approximately 18.7 lb.)		

Dimensions (Unit: mm)

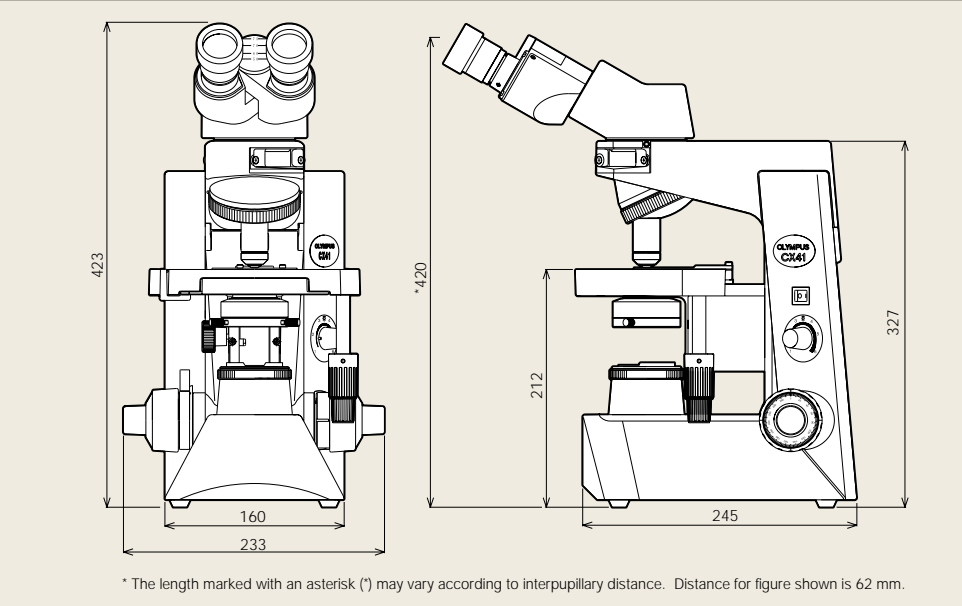

Objectives, Plan Achromat		
PLCN	Numerical Aperture (N.A.)	Working Distance (W.D.)
4X	0.10	18.5 mm
10X	0.25	10.6 mm
20X	0.4	1.2 mm
40X	0.65	0.6 mm
60X	0.8	0.2 mm
100XO	1.25	0.13 mm
100XOI	1.25–0.6	0.13 mm

Eyepiece	
	Field Number (F.N.)
WHB10X	20



ECO-PRODUCTS

CX41 is the environmental conscious product according to OLYMPUS' own standards.

Main features of OLYMPUS Eco-products are as follows.

- Lead-free and arsenic-free Eco-glass for optics, such as lenses and prisms.
- Exclusion of hexavalent chrome, mercury, lead and cadmium from metal materials and surface treatment of metal.
- Exclusion of lead solders.
- Adoption of cardboard for packing materials without styrene foam for promoting the recycling.

* A definition of exclusion depends on OLYMPUS standard. Some accessories are inapplicable.

Please visit our web site for further information: <http://www.olympus.co.jp/en/eco-products/>

- OLYMPUS CORPORATION is ISO14001 certified.
- OLYMPUS CORPORATION is FM553994/ISO9001 certified.
- Illumination devices for microscope have suggested lifetimes. Periodic inspections are required. Please visit our web site for details.
- Specifications and appearances are subject to change without any notice or obligation on the part of the manufacturer.



www.olympus.com

OLYMPUS CORPORATION
Shinjuku Monolith, 3-1, Nishi Shinjuku 2-chome, Shinjuku-ku, Tokyo, Japan
OLYMPUS EUROPA HOLDING GMBH
Wendenstrasse 14-18, 20097 Hamburg, Germany
OLYMPUS AMERICA INC.
3500 Corporate Parkway, Center Valley, Pennsylvania 18034-0610, U.S.A.
OLYMPUS SINGAPORE PTE LTD.
491B River Valley Road, #12-01/04 Valley Point Office Tower, Singapore 248373

OLYMPUS AUSTRALIA PTY. LTD.
31 Gilby Road, Mt. Waverley, VIC 3149, Melbourne, Australia.
OLYMPUS LATIN AMERICA, INC.
5301 Blue Lagoon Drive, Suite 290 Miami, FL 33126, U.S.A.
OLYMPUS (CHINA) CO., LTD.
A8F, Ping An International Financial Center, No. 1-3, Xinyuan South Road,
Chaoyang District, Beijing, China, 100027

OLYMPUS®

Your Vision, Our Future

System Microscope

CX41

CX2 Series


UIS2
World-leading optics

Optics and Performance Outstanding in its Class




Advanced optical and system performance with excellent cost-efficiency

The evolution of this bestselling microscope delivers new, cost-efficient improvements in both optics and system performance. Newly equipped with Olympus' leading-edge UIS2 optical system, it provides significantly enhanced image clarity in a variety of observation methods, from brightfield to reflected light fluorescence. For all inspection and training applications in the fields of biology and medicine, the CX41 sets the pace for its class in both basic and system performance.

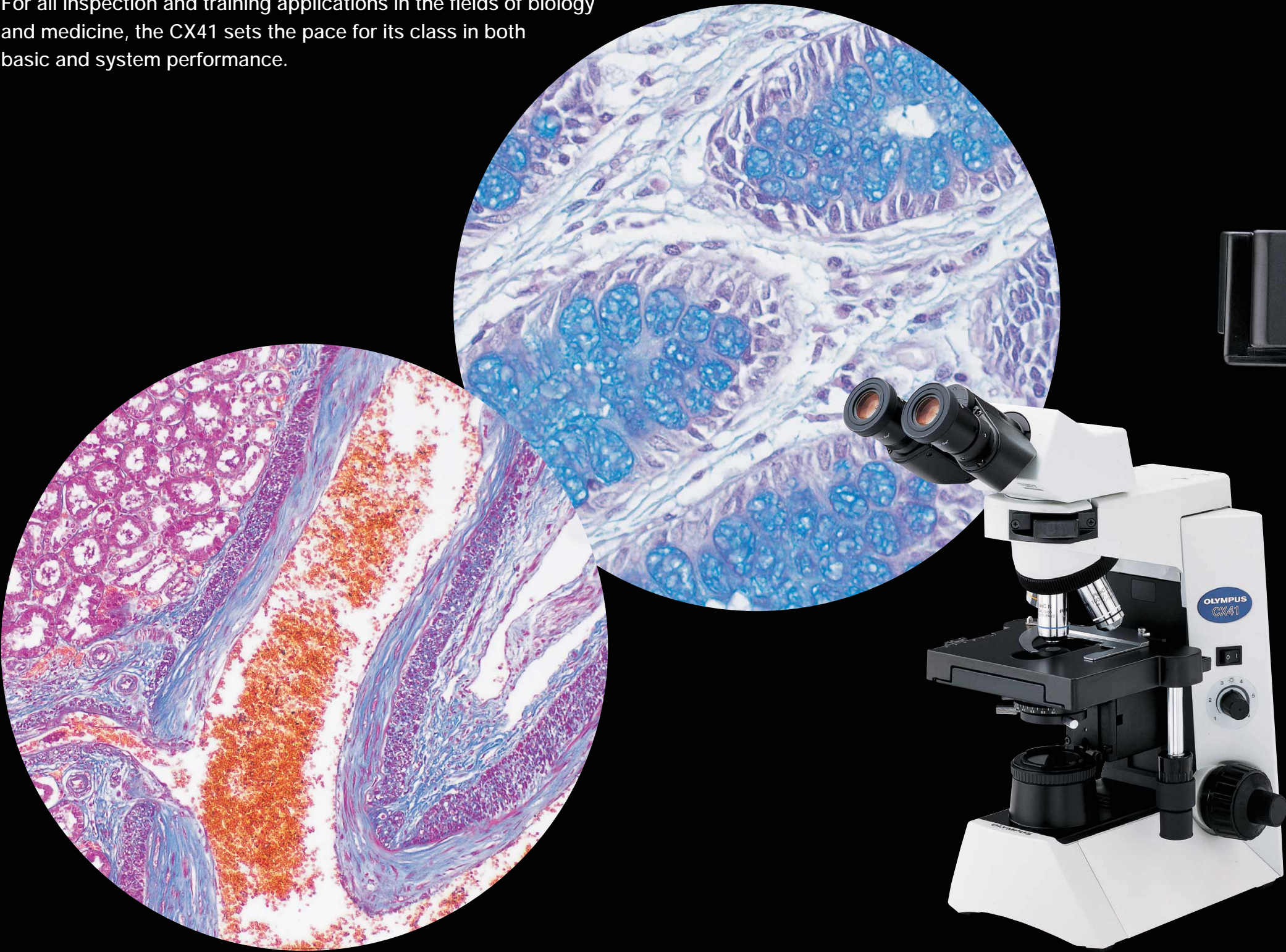

Outstanding flat images from PLCN objectives

The CX41 provides images of outstanding brightness and clarity in a variety of observation modes. As well as Olympus' renowned UIS2 optics infinity system, it employs the PLCN series of Plan Achromat objectives, which are made from carefully selected top quality glass and manufactured with the most rigorous precision. The result is a major improvement in image flatness, with the 10X and 40X objectives in particular providing images that are among the very best in this class of microscope. Transmitted light illumination is from a 6 V, 30 W high-intensity halogen light source.

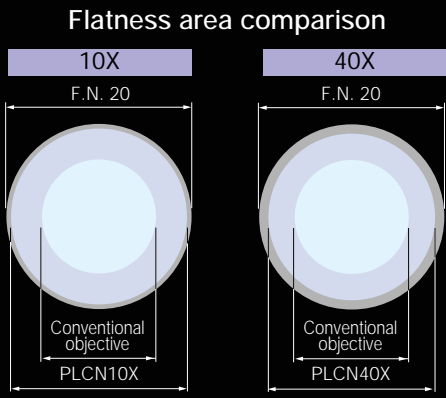

Excellent cost-performance in reflected light fluorescence and other observation methods



**Slide condenser / CX-SLC
Brightfield condenser / CH3-CD**
These Abbe type condensers allow brightfield observations from 4X to 100X. Accurate centering is provided by the attachment lens (CX-AL) and the iris diaphragm, to exclude unnecessary light and obtain bright Koehler illumination right across the magnification range. These highly economical condensers enable phase contrast and darkfield observations by simply adding basic accessories.



Simple phase contrast attachment / CX-PH1, 2, 3
For phase contrast observations at 10X, 40X and 100X.




Darkfield central stop / CH2-DS
For darkfield observations from 4X to 40X.
* Separate filter holder (CH2-FH) or attachment lens (CX-AL) required.



Low magnification adapter /CX-LA
Use of the 2X low magnification objective allows macro observation.

Reflected light fluorescence attachment / CX-RFA-2
Users can choose between blue or green excitation and transmitted light observations. UIS2 optics provide bright fluorescence images, with no intermediate magnifications when changing from transmitted light to fluorescence observation. Standard PLCN objectives can be used without replacement.

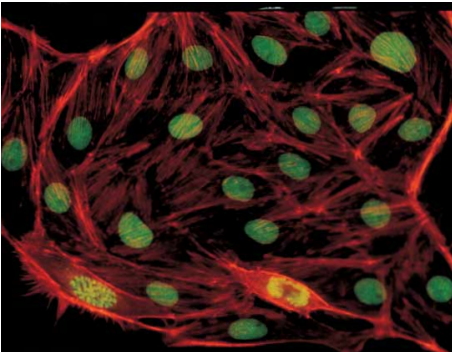

Simple polarizing condenser / CH3-CDP
With the optional plate adapter U-TAD, polarizing observations from 4X to 100X using a tint plate can be performed. A U-GAN analyzer is provided for gout inspection. Polarizing objectives from 4X to 100X are available.
* Separate polarizer U-POT and analyzer U-ANT required.



Dry darkfield condenser / CX-DCD
This dry-type darkfield condenser gives a superior darkfield effect without the need for immersion in oil. Suitable for use at 10X and 40X magnifications.



Phase-contrast condenser / CX-PCD
The multi-purpose CX-PCD condenser allows observation of brightfield, phase-contrast and darkfield images without exchanging condensers. Phase-contrast observation from 10x to 100x and darkfield observation from 10x to 40x is allowed.



Anti-fungus treatment
The treatment applied to the observation tubes, eyepieces and objectives, protects quality of optical parts even in high humidity regions.

Dependable basic performance assures outstanding operational convenience

Move the specimen with just one finger
Rubber grips are provided for the stage handles, allowing the specimen to be moved smoothly with just one finger. The slim body and conveniently positioned controls ensure that everything is within easy reach, so operators can maintain a natural posture.



Tilting binocular tube / U-CTBI
The tilting binocular tube lets each operator select the most suitable and comfortable eyepoint — a valuable contribution to reducing fatigue in extended observation sessions.



Inward-facing quintuple revolving nosepiece
The quintuple revolving nosepiece enables a wide range of magnification observations and various combinations of objectives. Unrestricted access to the space in front of the objective allows specimens to be exchanged quickly and easily.



Rackless stage with enhanced operability
To keep the work area clear, and to avoid interference with observation operations, the X-direction travel guide does not extend out from the side of the stage. The main and sub-scale displays are designed for easy read-out.



Torque adjustable focusing knob
The torque of the coarse focusing knob can be adjusted, to suit different operators' needs and to make focusing smooth and easy while keeping the hands on the desk. A stage upper limit stopper is also provided.



Easy transportation and installation
The CX41 is eminently portable, with convenient handgrips at the front and back of the frame and no inconvenient protrusion of the stage guide.



More accessories, more observation versatility



Dual observation attachment/ U-DO3
Enables dual, simultaneous observation of a single specimen from the same direction with equal magnification and brightness for both operators. A pointer can be used to indicate specific sections of the specimen to simplify the training process and enhance discussion.

Trinocular tube/ U-CTR30-2
Digital or video cameras can be attached for on-site/ remote illustration of educational/ discussion meetings.

Digital camera/ DP21
High-definition images displayed with stunning detail in real time. Seamless conference support with real-time image display and smooth operation.



CX41+DP21

Eyepoint adjuster/ U-EPA2
Allows the eyepoint position to be raised by 30 mm. Up to two eyepoint adjusters can be attached between the frame arm and observation tube.



Arrow pointer/ U-APT
Enables insertion of an LED arrow for display in a digital image.



2x magnification changer/ U-ECA
Magnification is doubled by engaging the auxiliary 2x lens.



Drawing attachment/ U-DA
The drawing attachment projects an image of the pencil and drawing surface into the visual field.



Trinocular intermediate attachment/ U-TRU
A binocular tube on its own allows digital imaging in combination with this trinocular intermediate attachment.

